



# PROSTATE CANCER BIOREPOSITORY NETWORK

SOP No: 005  
QC of DNA

---

STANDARD OPERATING PROCEDURE	<b>SOP No. 005</b> <b>QC of DNA</b>
<b><i>UNAUTHORIZED COPYING PROHIBITED</i></b>	<b>Version Number: 2.0 (November 2016)</b> Number of Pages: 3
Electronic Filename: PCBN.SOP05.v2.0 QC of DNA	



# PROSTATE CANCER BIOREPOSITORY NETWORK

SOP No: 005  
QC of DNA

## 1. PURPOSE

To describe the procedure for the QC of DNA extracted from frozen tissue by real-time PCR. NOTE: This SOP does not cover detailed procedures for handling Human Biological Materials or hazardous chemicals and it is recommended that personnel following this SOP refer to institutional safety guidelines.

**The following SOP for QC of DNA is followed by the JHU, NYU, MSKCC and UW Network Sites of the PCBN.**

## 2. RESPONSIBILITIES

**Authorized personnel** performing QC of DNA extracted from frozen prostate tissue must ensure that:

- all procedures are followed correctly
- all samples are adequately coded during processing
- all documentation is completed, and accurate records maintained on all samples

## 3. HEALTH AND SAFETY

Personnel carrying out this procedure must maintain safe working practices and observe all relevant Environmental Health and Safety (EH&S) guidelines. This includes the appropriate use of Personal Protective Equipment (PPE), and the procedures for waste disposal, disinfection & spill clean-up and biosafety.

## 4. EQUIPMENT AND MATERIALS FOR JHU & NYU NETWORK SITES

Equipment	Materials
iCycler real time Thermocycler Nanodrop	96-well plates Primers as described SYBR Green PCR Master Mix (Invitrogen Catalog # 4309155)

## 5. EQUIPMENT AND MATERIALS FOR UW & MSKCC NETWORK SITES

Equipment	Materials
Gel electrophoresis system Gel Doc Nanodrop	Agarose TAE Buffer 6X Sample Loading Buffer DNA Ladder Standard



---

## 6. PROCEDURES FOR JHU AND NYU NETWORK SITES ONLY

- Quantitate DNA samples extracted using **PCBN SOP004 DNA Extraction from Frozen Tissue** by Nanodrop. Refer to the Nanodrop user guide for instructions of use.
- If DNA concentration is greater than 100 ng/ul, dilute DNA sample to 100 ng/ul.
- QC of DNA samples extracted using **PCBN SOP004 DNA Extraction from Frozen Tissue** by real-time PCR.

- Rxn mix (20 µl): µl/rxn

gDNA	2
SYBR Green	10
F-primer	1
R-primer	1
dH <sub>2</sub> O	6

- A. β-globin primers (100 bp product)

Forward: 5'-GTGCACCTGACTCCTGAGGAGA-3'  
Reverse: 5'-CCTTGATACCAACCTGCCAG-3'

- B. 18S primers (110 bp product)

Forward: 5'-GATGGTAGTCGCCGTGCC-3'  
Reverse: 5'-GCCTGCTGCCTTCCTTGG-3'

- Cycling parameters

Cycle 1: ( 1X)  
Step 1: 95.0°C for 03:00  
Cycle 2: ( 35X)  
Step 1: 95.0°C for 00:30  
Step 2: 60.0°C for 00:30  
Step 3: 72.0°C for 00:45

## 7. PROCEDURES FOR UW AND MSKCC NETWORK SITES ONLY

- Quantitate DNA samples extracted using **PCBN SOP004 DNA Extraction from Frozen Tissue** by Nanodrop. Refer to the Nanodrop user guide for instructions of use.
- Measure the  $A_{260}/A_{280}$  ratio for an estimate of DNA purity with respect to contaminants that absorb UV.
- DNA length can be determined using agarose gel electrophoresis.
- DNA samples with sharp 10kb band are considered as high-quality DNA.